

Quick-Guide for Software Ver. 3.0

Revision 002 – 13th July 2016 – Build on Systemvars 94.102

Gram Commercial A/S
Aage Grams Vej
6500 Vojens, Danmark
Tel: +45 73 20 13 00
Fax: +45 73 20 12 01
www.gram-bioline.com


Biostorage you can depend on

Quick guide for software version 3.0

Cabinet On/OFF

Press the Q-key to switch on the cabinet. Press and hold it for 6 sec. the cabinet will switch off (Stand-by). During the start-up the program variant and SW version will be presented shortly. Revision number can be presented under "Display test".

Other short cuts

Using the key-pad functions is accessed or activated.

Key combination:	Time:	Function:
P+Q	> 3 sec.	Start or stop a defrosting cycle manually.
Q+1	> 6 sec.	Key pad lock on/off. With a lock key it's not possible to access any menu or make any changes.
(+)	-	Showing highest recorded temperature peak since the last reset of alarm history.
(-)	-	Showing lowest recorded temperature peak since the last reset of alarm history.
(+) + (-)	> 6 sec.	Re-setting alarm history.
P+1+3	> 6 sec.	Re-setting all changes in the parameters and brings the settings back to default.
P+1	> 6 sec.	Access to User menu and alarm settings . See more in the tables below.
P+2	> 6 sec.	Access to System monitoring and presentations . See more in the tables below.
P+3	> 6 sec.	Access to Cooling system setup . See more in the tables below.
P+4	> 6 sec.	Access to the Test Program . See more in the tables below.
P+5	> 6 sec.	Access to the Control of the Sensors . See more in the tables below.

Changes in product setup

If changes are needed in the controllers settings, then enter the menus as according to above. The values is changed using the + and - keys. The new setting is saved by pressing the P-key. Leave the menus with Q-key.

User menu and alarm settings

The alarm system is separated into two. One alarm system triggers alarms only locally, which means the error only is displayed. The second alarm system triggers locally and external. Error codes in the display, but also alarm activates also the potential-free alarm output. Each alarm system is working independently.

Menu access P+1 →	↓	→		
Dry cooling	dC			Activation of dry cooling program. [H1=off / H0=active]
Local alarm setting	LAL	LhL	[° C]	Setting the upper alarm limit. At alarm the display is showing: [A2].
		LLL	[° C]	Setting the lower alarm limit. At alarm the display is showing: [A3].
		Lhd	[min.]	Time delay for upper alarm limit.
		LLd	[min.]	Time delay for lower alarm limit.
		dA	On/off	Activation of door alarm [1=on / 0=off]. At alarm the display is showing: [A1].
		dAd	[min.]	Time delay for door alarm.
		BU	On/off	Activation of buzzer [1=on / 0=off]. The buzzer sounds at alarms [A1], [A2], [A3].
External alarm setting	EAL	EhL	[° C]	Setting the upper alarm limit. At alarms the display is showing: [A4].
		ELL	[° C]	Setting the lower alarm limit. At alarms the display is showing: [A5].
		Ehd	[min.]	Time delay for upper alarm limit.
		ELd	[min.]	Time delay for lower alarm limit.
		dA	On/off	Activation of door alarm [1=on / 0=off]. At alarms the display is showing: [A1].
		dAd	[min.]	Time delay for door alarm.
		BU	On/off	Activation of buzzer [1=on / 0=off]. The buzzer sounds at alarms [A1], [A4], [A5].
Sensor calibration	cAL	cA	[° K]	Offset adjustment on A-sensor. Cabinet sensor.
		cE	[° K]	Offset adjustment on E-sensor. Extra sensor (placed in air or in bottle).
		cF	[° K]	Offset adjustment on F-sensor. Sensor for frost protection.
Frost protection	FP	Act	On/off	Activation on frost protection.
		tES	On	Test the frost protection. Switch off compressor after time = C4.
		SEt	[° C]	Adjustment of cut-out for compressor.
		PrE	[...]	Presentation of F-sensor.
	ALL			Activation of escorting alarm limits. [FAS] = locked borders / [ESC] = following set point.
	dEF			Numbers of defrosting pr. day.
	dPS			Selection of sensor displaying in the display. Choose between: A, E or F.
The alarms:				Text in display and description:
Alarm messages from above.	A1			Door alarm is triggered either by the LAL or EAL alarm systems.
	A2			Alarm code triggered by a too hot temperature associated with the local alarm system.
	A3			Alarm code triggered by a too cold temperature associated with the local alarm system.
	A4			Alarm code triggered by a too hot temperature associated with the external alarm system.
	A5			Alarm code triggered by a too cold temperature associated with the external alarm system.
	A6			Frost protection has stopped compressor and ensured that the room temperature is not too cold.

Temperature monitor and presentations

Menu for setting the monitoring, alarm history, sensor choice for alarm systems and temperature read-out.

Menu access P+2 →	↓	→		
Alarm setting	A	A1	[° C]	By overheated condenser this starts the compressor protection program
		A2	[° C]	Temperature in condenser to de-activate the compressor protection program.
		A4	On/off	Activation of alarm history [1=on / 0=off]. The program only works with alarm system: LAL .
		A5		Choose between A, E or F sensor for the alarm systems LAL and EAL .
Presentation in display	P	P1	[min.]	Freezing the temperature in minutes in display after a defrosting cycle.
		P2	[° K]	Freezing the temperature in display under normal run. Temperature fluctuation filter.
		P3	[sec.]	Display refresh rate.
		P4		Selection between Celsius or Fahrenheit temperature scale.

Settings for cooling system

Menu for setting the evaporator fan, compressor and defrosting cycles.

Menu access P+3 →	↓	→	RX-model V1+	BioUltra E3+	
Compressor setting	C	C1		C1	[° K] Differential for compressor cut-in and cut-out.
		C2		C2	[° C] Maximum allowable temperature limit.
		C3		C3	[° C] Lowest allowable temperature limit.
		C4		C4	[min] Forced break time for the compressor between cut-out and cut-in
		C5			Number of condenser sensor connected to controller
		C6			[min] Permitted time the door can be open before the compressor stops.
			C7		[° K] (RX-model / Soft differential for cool/heat cut-out).
				C8	[° C] (BioUltra / Set point for condenser fan.)
				C9	[° K] (BioUltra / Cut-out differential for condenser fan).
Evaporator fan setting	F	F1		L1	[° C] Temperature for fan start after the defrosting cycle. Start of LT compressor in E3+
		F2			[min.] Time evaporator is paused while the compressor is stopped.
		F3			[sec.] Time evaporator fan is in operation while the compressor is stopped
				L4	[° C] Stop temperature in cascade heat exchanger for LT compressor – At normal run.
				L5	[° C] Stop temperature in cascade heat exchanger for LT compressor – By abnormal run.
				L6	[° K] Temperature deviation for choosing between "normal or abnormal" run.
				L7	Running mode for LT compressor [1=slave / 0=independent].
Defrosting cycle setting	d	d1			Numbers of defrosting pr. day.
		d2			[° C] Defrosting stop temperature measured in the evaporator
		d3			On/off Activating the defrosting cycle on first power-up start [0=on / 1=off]
		d4			[min.] Maximum defrosting time.
		d5			Choose between automatic [1], air [2] or electrical defrosting method [3]
		d6			[min.] Dripping time after defrosting cycle ends.
		d7			[° C] The maximum level that determines the defrosting method when d5 = [1]
		d8			[° C] Temperature in evaporator to start an extra defrosting cycle automatically.
				d10	[min.] Defrosting window.
				d11	[h] Time delay at 100 % compressor running before forced defrosting cycle.

Test program

Individually control of relay outputs and connected external components. Select the menu item and press the P-key. The relay switches and power is connected to external component. The display shows [on]. Stop the test with Q-key.

Menu access P+4 →	↓	→	BioUltra E3+	P-key → [on] / Q-key [off]
Compressor	tC	tHt		Control of compressor and condenser fan / (BioUltra: Control of HT compressor).
Evaporator fan	tF	tLt		Control of evaporator fan / (BioUltra: Control of LT compressor).
Defr. heating element	td	td		Control of defrosting heating element. Warning: The heating element becomes very hot. Danger!
Light	tL	tCF		Control of light / (BioUltra: Control of condenser fan speed change – 900 rpm to 1200 rpm).
Alarm output	tA	tA		Control of the potential free relay out-put.
Display test	tdP	tdP		All LED's in the display will lights up for 1 sec. followed by the software revision number.

Control of sensor and display

With this menu it's possible to have instantaneous sensor read-outs.

Menu access P+5 →	↓	P-key → [° C]	Message in display and cause
Cabinet sensor	P-A		F1 Error on cabinet sensor
Evaporator sensor	P-b		F2 Error on evaporator sensor
Condenser sensor 1	P-C		F3 Error on condenser sensor 1
Condenser sensor 2	P-d		F4 Error on condenser sensor 2
Extra/Reference sensor	P-E		F5 Error on Extra/Reference sensor
Sensor for frost protect.	P-F		F6 Error on sensor for frost protection
Overheated condenser caused by a clogged grease filter. Triggered by both C and d sensor			
BioUltra: High pressure switch has triggered by excessive pressure in the LT cooling system.			
When the door opens the symbol lights up. A too long open door this will trigger the alarm [A1].			
			F7 Overheated condenser 1 or 2.
			F8
			-0- Symbol for open door



Biostorage you can depend on

Default factory settings

Below are all BioLine program variants shown. In the columns under each program variant, the factory settings is presented. Should any questions occur, then please take contact to BioLine Tech Support.

		Refrigeration						Extended		Freezer				Heat		Extreme low temp. range				
Systemware revision number		K2+	K4+	K5+	K6+	K8+	K9+	K70	M4+	M5+	F5+	F6+	F51	F61	F70	V1+	E1+	E3+	E4+	
		94.102	94.102	94.102	94.102	94.102	94.102	94.98	94.80	94.102	94.80	94.80	94.80	94.80	94.98	94.80	94.80	94.80	94.80	
Setpoint (celsius)		5	5	5	5	5	5	5	4	5	-20	-20	-20	-20	-20	5	-35	-80	-40	
Temperature range (celsius)		+20/+2	+20/+2	+20/+2	+20/+2	+20/+2	+20/+2	+15/+2	+6/+2	+20/+2	-5/-25	-5/-25	-5/-25	-5/-25	-5/-25	+45/0	-5/-38	-60/-95	-5/-45	
User menu: P=1																				
Dry refrigeration on=H0/Off=H1		dC																		
Thawing speed		UF																		
Local Alarm Limits		LAL																		
Upper local alarm limit (celsius)		UHL						ELH		F5H				F70H		V1H				
Lower local alarm limit (celsius)		LLH						ELL		F5L				F70L		V1L				
Time delay for upper alarm limit (min.)		LHd						ELd		F5d				F70d		V1d				
Time delay for lower alarm limit (min.)		LLd						ELd		F5d				F70d		V1d				
Door alarm on=1/Off=0		dA						dA		dA				dA		dA				
Time delay for open door (min.)		dAd						dAd		dAd				dAd		dAd				
Buzzer for local alarm on=1/Off=0		bU						bU		bU				bU		bU				
External Alarm Limits		EAL																		
Upper external alarm limit (celsius)		EHL						EHL		E5H				E70H		V1H				
Lower external alarm limit (celsius)		ELL						ELL		E5L				E70L		V1L				
Time delay for upper alarm limit (min.)		EHLd						EHLd		E5d				E70d		V1d				
Time delay for lower alarm limit (min.)		ELld						ELld		E5d				E70d		V1d				
Door alarm on=1/Off=0		dA						dA		dA				dA		dA				
Time delay for open door (min.)		dAd						dAd		dAd				dAd		dAd				
Buzzer for local alarm on=1/Off=0		bU						bU		bU				bU		bU				
Calibration of sensor		CAL																		
Offset adjustment sensor A (kwh)		CA						CA		CA				CA		CA				
Offset adjustment sensor F (kwh)		CF						CF		CF				CF		CF				
Offset adjustment sensor E (kwh)		CE						CE		CE				CE		CE				
Frost Protection		FP																		
Activation of frost protection On=1/Off=0		Act																		
Test of frost protection		IES						IES		IES				IES		IES				
Setpoint of frost protection (celsius)		SEt						SEt		SEt				SEt		SEt				
Display of current sensor temperature (celsius)		PRE						PRE		PRE				PRE		PRE				
Fixed or escaping alarm limits (FAS - ESC)		ALL						SCL		FAS				FAS		FAS				
Soft chilling (soft-chl)		SCL						SCL		SCL				SCL		SCL				
Hard chilling (hard-chl)		HCL						HCL		HCL				HCL		HCL				
Time controlled chilling (timed-chl)		PCL						PCL		PCL				PCL		PCL				
Number of defrosts / 24h		dF						dF		dF				dF		dF				
Selected sensor display		dPS						dPS		dPS				dPS		dPS				
Possible choice for sensor displaying in display		A,E,F						A,E,F		A,E				A,E		A,E				
Alarm & presentation: P=2		A																		
Alarm settings		A																		
Condenser monitoring alarm on (celsius)		A1						A1		A1				A1		A1				
Condenser monitoring alarm off (celsius)		A2						A2		A2				A2		A2				
Geninndrædes liden for akustisk alarm (min.)		A3						A3		A3				A3		A3				
Alarm history on=1/Off=0		A4						A4		A4				A4		A4				
Selection of sensor for alarm system		AA						AA		AA				AA		AA				
Possible choice for sensor displaying in display		A,E,F						A,E,F		A,E				A,E		A,E				
Presentation of temperature		P																		
Temperature display hold after defrost		P1						P1		P1				P1		P1				
Temperature display hold during normal operation		P2						P2		P2				P2		P2				
Display updating frequency (sec)		P3						P3		P3				P3		P3				
Temperature display: Celsius or Fahrenheit		P4						P4		P4				P4		P4				
System setup: P=3		S																		
Compressor settings		C																		
Differential for compressor start and stop (kwh)		C1						C1		C1				C1		C1				
Max. allowed setpoint (celsius)		C2						C2		C2				C2		C2				
Min. allowed setpoint (celsius)		C3						C3		C3				C3		C3				
Compulsory compressor pause time (min.)		C4						C4		C4				C4		C4				
No. of sensors for condenser monitoring		C5						C5		C5				C5		C5				
Door open time before compressor stops (min.)		C6						C6		C6				C6		C6				
Soft differential for cool/heats cut-out (kwh)		C7						C7		C7				C7		C7				
Setpoint for condenser fan		C8						C8		C8				C8		C8				
Cut-out differential for condenser fan (kwh)		C9						C9		C9				C9		C9				
Evaporator fan settings		F																		
Evap. fan start after defrost, and during dry ref. (celsius)		F1 (L1)						F1		F1				F1		F1				
Evap. fan pause time at compressor stop (min.)		F2						F2		F2				F2		F2				
Ford. vent. start ved kompressor stop (sek.)		F3						F3		F3				F3		F3				
Stop temp. of LT compressor (in cascade evaporator at normal deviation (celsius)		F4 (L4)						F4		F4				F4		F4				
Stop temp. of LT compressor in cascade evaporator at large deviation (celsius)		F5 (L5)						F5		F5				F5		F5				
Temp. deviation for determining between "normal" or "abnormal" run		F6 (L6)						F6		F6				F6		F6				
Operation mode for LT compressor (slave or independent)		F7 (L7)						F7		F7				F7		F7				
Defrost settings		d																		
Number of defrosts / 24h		d1						d1		d1				d1		d1				
Stop temperature in evaporator during defrost (celsius)		d2						d2		d2				d2		d2				
Defrosting on/off = (p1) at power up first time		d3						d3		d3				d3		d3				
Max allowed defrosting time (min.)		d4						d4		d4				d4		d4				
Defrosting mode (1=automatic, 2=air, 3=electrical)		d5						d5		d5				d5		d5				
Dip time after defrost stop (min.)		d6						d6		d6				d6		d6				
Limit for automatic defrosting mode (celsius)		d7						d7		d7				d7		d7				
Evaporator monitoring (celsius)		d8						d8		d8				d8		d8				
Defrost after chilling stop		d9						d9		d9				d9		d9				
Defrosting window after compressor cut-out (min.)		d10						d10		d10				d10		d10				
Time for overruled defrosting cycle at 100 % compressor running (hours)		d11						d11		d11				d11		d11				
Sensor configuration		St																		
St. = +25 °C / 1505 = 40 °C / 1513 = -80 °C		St																		
St. = +25 °C / 1505 = 40 °C / 1513 = -80 °C		St																		
St. = +25 °C / 1505 = 40 °C / 1513 = -80 °C		St																		
St. = +25 °C / 1505 = 40 °C / 1513 = -80 °C		St																		
St. = +25 °C / 1505 = 40 °C / 1513 = -80 °C		St																		
Other Software Features		dC																		
Dry Cooling		UF																		
Thaw-out		UF																		
Blaschiller		UF																		
Compressor protection program		UF																		
Demand controlled condenser fan controlled		UF																		
Emergency program at defective room sensor		UF																		
Evaporator surveillance program		UF																		
Alarm system designed for Commercial use		UF																		
Alarm system designed for BIOBASIC		UF																		
Simple alarm system designed BioCompact		UF																		
Advanced alarm system designed for BioLine & Biocompact II		UF																		
Alarm history feature		UF																		
Low temperature protection connectivity		UF																		
Presenting sensor A and E in display		UF																		
Presenting sensor A and F in display		UF																		
Presenting sensor A, E and F in display		UF																		
Current display of temperature		UF																		
No evaporator sensor viable		UF																		
Defrosting dot in display at defrosting cycle		UF																		
Manual On/Off operation for light in glass door		UF																		
Automatically cut-out for traditional light after 30 minutes		UF																		
Indicator activated at low temperature protection alarm		UF																		
ASDD - Delayed defrosting cycle depending on running compressor		UF																		
Automatically Start-up at power-On		UF																		





LABORATORY EQUIPMENT MAINTENANCE, REPAIR, CALIBRATION AND SALES

Established in 1987, Henderson Biomedical is the UK's leading laboratory equipment sales and service provider. Our knowledgeable team can provide you with excellent sales advice on a range of different types of laboratory equipment including centrifuges, refrigerators, freezers and heat sealers.

Henderson Biomedical is also able to provide you with first class after-sales service and calibration of your laboratory equipment. We are an **ISO 17025 (UKAS) accredited calibration test laboratory** and our team of Field Service Engineers cover the whole of the United Kingdom.

Please contact us for more information on the types of equipment we supply and the different after-sales services we can offer.

Henderson Biomedical

Unit 3, Swan Close
Croydon CR0 2DZ
United Kingdom

Tel: 020 8663 4610

For sales enquiries: sales@henderson-biomedical.co.uk
For all other enquiries: info@henderson-biomedical.co.uk
www.henderson-biomedical.co.uk

