

# ROTOFIX 32 A

## Rugged and indispensable

For decades, the ROTOFIX 32 A has set the standard in daily lab routine thanks to its versatility and solid construction. The benchtop centrifuge spins sample volumes up to 4 x 100 ml, 40 blood collection tubes or 8 x 50 ml conical tubes with a simple user interface. Hettich's cytology rotors are compatible with most existing funnel/slide systems and have bio-containment lids.

### — FEATURES

- RPM: 500 - 6,000 min<sup>-1</sup> – adjustable in increments of 100
- Max. RCF: 4,226
- Max. capacity: 4 x 100 ml / 6 x 85 ml
- Choice of 9 rotors
- IVDR-conform according to regulation (EU) 2017/746
- Easy operation with keypad
- 2 individual deceleration stages

### — FIELDS OF APPLICATION

- Hospitals
- Cell culture laboratories
- University / Academic research
- Pharmaceutical laboratories
- Physician's Office Laboratory (POL)



Centrifuge packages for the model can be found on [page 65](#)



More information about the control panel can be found on [page 202](#)

**CYTO**

Cyto system available for this model. More information on [page 180](#)



according to regulation (EU) 2017/746

## TECHNICAL DATA

<b>ROTOFIX 32 A</b>	
voltage *)	208 – 240 V 1 ~
frequency	50 – 60 Hz
consumption	300 VA
emission, immunity	EN/IEC 61326-1, class B
max. capacity	4 x 100 ml / 6 x 94 ml
max. RPM	6,000 min <sup>-1</sup>
max. RCF	4,226
running time	1–99 min, ∞ continuous run, short cycle mode (impulse button)
dimensions (WxDxH)	366x430x257 mm
weight	approx. 23 kg
noise level	52 dB (A) with rotor 1628
<b>Cat. No.</b>	<b>1206</b>

100 – 127 V 1 ~ / 50 – 60 Hz \*)

1206-01







emission, immunity

FCC class B




\*) Other voltages on request.
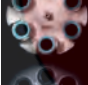
## AVAILABLE ROTORS

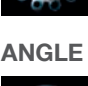
### SWING-OUT ROTORS

	angle	max. RPM	max. capacity	Cat. No.	page
 swing-out rotor, 4-place	90°	4,000 min <sup>-1</sup>	4x100 ml	<b>1624</b>	56
 swing-out rotor, 4-place	90°	4,000 min <sup>-1</sup>	4x100 ml	<b>1324</b>	59
 swing-out rotor, 6-place	90°	4,000 min <sup>-1</sup>	6x50 ml	<b>1619</b>	61
 swing-out rotor, 8-place	45°	4,000 min <sup>-1</sup>	8x50 ml	<b>1617</b>	61
 swing-out rotor, 8-place	90°	4,000 min <sup>-1</sup>	8x15 ml	<b>1611</b>	63
 swing-out rotor, 12-place	55° / 60° / 80°	4,000 min <sup>-1</sup>	12x15 ml	<b>1628</b>	63

### ANGLE ROTORS

 angle rotor, 8-place	45°	4,000 min <sup>-1</sup>	8x50 ml	<b>1418</b>	63
 angle rotor, 6-place	35°	6,000 min <sup>-1</sup>	6x85 ml	<b>1620A</b>	65
 angle rotor, 12-place	35°	6,000 min <sup>-1</sup>	12x15 ml	<b>1613</b>	65

## SWING-OUT ROTOR, 4-PLACE | 1624

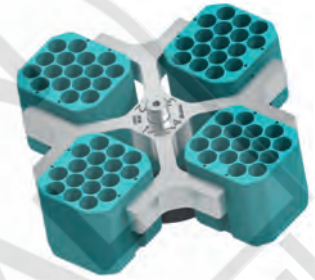

### Rotor

max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,451
max. capacity	4 x 50 ml
run-up   run-down, braked in sec	22   25
angle   max. noise level	90°   55 dB (A)
<b>Cat. No.</b>	<b>1624</b>

### Vessels

capacity in ml	5	5	6	7	9	9	15	15	20	25	45	50
Ø x L in mm	12x75	12x75	12 x 82	12 x 100	14 x 100	14 x 100	17 x 100	17 x 100	21 x 100	24 x 100	31 x 100	34 x 100
max. RCF <sup>2)</sup>	2,218	2,164	2,343	2,343	2,308	2,415	2,308	2,451	2,361	2,451	2,361	2,451
radius in mm	124	121	131	131	129	135	129	137	132	137	132	137

<b>Cat. No.</b>	<b>tubes<sup>3)</sup></b>											
<b>+</b>	with decanting aid		with decanting aid	with decanting aid		+ 0701						
<b>Carrier</b>												
boring Ø x L in mm	12 x 75	13.5x65	12.5x71.5	12.5x71.5	14.6x74	14.6x78	17.6x71.5	17.6x78	21.5x74	26x78	32x74	35x78
vessels per rotor	16	68	16	16	20	40	16	28	8	8	4	4
<b>Cat. No.</b>	<b>1369-91</b>	<b>1372</b>	<b>1369-92</b>	<b>1369-92</b>	<b>1370</b>	<b>1741</b>	<b>1369</b>	<b>1742</b>	<b>1346</b>	<b>1745</b>	<b>1345</b>	<b>1746</b>

### Vessels

capacity in ml	1.1-1.4	2.6-3.4	4-5.5	4.9	1.6-5	4-7	4-7	8.5-10	30	1-8
Ø x L in mm	8 x 66	13 x 65	15 x 75	13 x 90	13 x 75	16 x 75	13 x 100	16 x 100	26 x 95	simple / multiple
max. RCF <sup>2)</sup>	2,415	2,325	2,325	2,451	2,325	2,325	2,451	2,451	2,451	1,646
radius in mm	135	130	130	137	130	130	137	129	137	92

<b>Cat. No.</b>	<b>blood collection / urine vessels</b>								-	<b>cyto chambers</b>
<b>+</b>	+ 0701	+ 0716	+ 0716		+ 0716	+ 0716				
<b>Carrier</b>										
boring Ø x L in mm	14.6 x 78	17.6 x 78	17.6 x 78	14.6 x 78	17.6 x 78	17.6 x 78	13.5 x 78	17.6 x 71.5	26 x 78	-
vessels per rotor	40	28	28	40	28	28	28	16	8	4
<b>Cat. No.</b>	<b>1741</b>	<b>1742</b>	<b>1742</b>	<b>1741</b>	<b>1742</b>	<b>1742</b>	<b>1739</b>	<b>1369<sup>4)</sup></b>	<b>1745</b>	<b>1660</b>

**CYTO** Cyto system available for this model. More information on page 180

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.  
 4) Please remove the spacer.  
 16) Packed in units of 10 pieces.  
 20) Vacutainers made of glass may not be used.

## SWING-OUT ROTOR, 4-PLACE | 1624



### Rotor

max. RPM   max. RCF	4,000 min <sup>-1</sup>   1,968
max. capacity	48 x 4 ml
run-up   run-down, braked in sec	22   25
angle   max. noise level	90°   56 dB (A)
<b>Cat. No.</b>	<b>1624</b>

### Bucket

Lid bioseal <sup>5)</sup>	1492
<b>Cat. No.</b>	<b>1366</b>



### Vessels

capacity in ml	1.5	2.0	1	3	4
Ø x L in mm	11 x 38	11 x 38	6 x 45	10 x 60	12 x 60
max. RCF <sup>2)</sup>	1,968	1,968	1,950	1,932	1,932
radius in mm	110	110	109	108	108
<b>Cat. No.</b>	<b>microliter tubes</b>	<b>Rhesus</b>	<b>tubes<sup>2)</sup></b>		



### Adapter

boring Ø x L in mm	11.5 x 38	11.5 x 38	6.5 x 23	10.5 x 23	12.5 x 44
vessels per rotor	36	36	120	48	48
<b>Cat. No.</b>	<b>5277</b>	<b>5277</b>	<b>1357</b>	<b>1327</b>	<b>1326</b>



## SWING-OUT ROTOR, 4-PLACE | 1624




### Rotor

max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,665
max. capacity	4 x 100 ml
run-up   run-down, braked in sec	22   25
angle   max. noise level	90°   55 dB (A)
<b>Cat. No.</b>	<b>1624</b>


### Bucket

Lid bioseal <sup>5)</sup>	1492
<b>Cat. No.</b>	<b>1481</b>

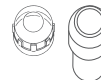


### Vessels

capacity in ml	0,5	1,5	2,0	1	3	5	6	7	9	15	25	50	85	100
Ø x L in mm	10,7 x 46	11 x 38	11 x 38	6 x 45	10 x 60	12 x 75	12 x 82	12 x 100	14 x 100	17 x 100	24 x 100	34 x 100	38 x 101	44 x 100
max. RCF <sup>2)</sup>	2,379	2,451	2,451	2,594	2,630	2,558	2,558	2,558	2,540	2,540	2,433	2,415	2,612	2,558
radius in mm	133	135	135	145	147	141	141	141	142	142	136	135	146	143
<b>Cat. No.</b>	<b>Pediatric</b>	<b>microtubes</b>		<b>Rhesus</b>	<b>tubes<sup>2)</sup></b>									

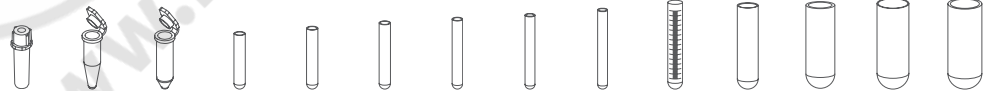

### Adapter

boring Ø x L in mm	11.2 x 38	11.2 x 38	11.2 x 38	6.5 x 34	10.5 x 43	13.4 x 48	13.4 x 48	13.4 x 48	17.6 x 89	17.6 x 89	25.2 x 77	35.2 x 77	38.5 x 92	45.9 x 98
vessels per rotor	20	20	20	108	36	20	20	20	16	16	4	4	4	4
<b>Cat. No.</b>	<b>1351</b>	<b>1351</b>	<b>1351</b>	<b>1339</b>	<b>1343</b>	<b>1383</b>	<b>1383</b>	<b>1383</b>	<b>1329</b>	<b>1329</b>	<b>1330</b>	<b>1331</b>	<b>1396</b>	<b>0761</b>

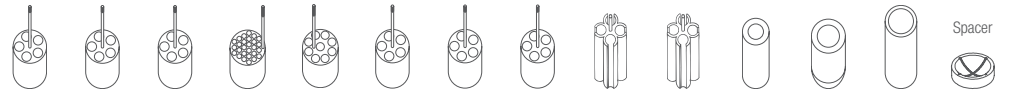

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.  
 5) Tested by TÜV in conformity with DIN EN 61010, section 2-020.



## SWING-OUT ROTOR, 4-PLACE | 1624




### Rotor

max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,665
max. capacity	4 x 100 ml
run-up   run-down, braked in sec	22   25
angle   max. noise level	90°   55 dB (A)
<b>Cat. No.</b>	<b>1624</b>

### Bucket

Lid bioseal <sup>5)</sup>	1492
<b>Cat. No.</b>	<b>1481</b>

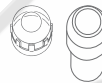

### Vessels

capacity in ml	1.1–1.4	2.7–3	2.6–4.9	4–8.5	9–10	10	1.6–7	4–10	15	50	12	25	30	50	10
Ø x L in mm	8 x 66	11 x 66	13 x 65/90	15 x 75/92	16 x 92	15 x 102	13 x 75/100	16 x 75/100	17 x 120	29 x 115	17 x 100	25 x 90	25 x 110	29 x 115	16 x 80
max. RCF <sup>2)</sup>	2,540	2,558	2,558	2,576	2,540	2,665	2,558	2,522	2,665	2,665	2,665	2,343	2,665	2,665	2,522
radius in mm	142	143	143	144	142	149	143	141	149	149	149	131	149	149	141
<b>Cat. No.</b>	<b>blood collection / urine vessels</b>								<b>tubes with screw cap</b>						



### Adapter

boring Ø x L in mm	9 x 49	13.4 x 48	13.4 x 48	15.6 x 47	17.6 x 91	17.6 x 91	13.4 x 48	16.5 x 56	17 x 90	30 x 90	17 x 80	26 x 72	26 x 80	29.5 x 80	16.5 x 56
vessels per rotor	28	20	20	16	16	16	20	16	4	4	4	4	4	4	16
<b>Cat. No.</b>	<b>1457</b>	<b>1383</b>	<b>1383</b>	<b>1459</b>	<b>1329</b>	<b>1329<sup>4)</sup></b>	<b>1383</b>	<b>1348</b>	<b>1347</b>	<b>1384</b>	<b>6311</b>	<b>1363</b>	<b>1365</b>	<b>6318</b>	<b>1348</b>

### Vessels

capacity in ml	30	50	85	30
Ø x L in mm	26 x 95	29 x 107	38 x 106	44 x 105
max. RCF <sup>2)</sup>	2,451	2,630	2,612	2,540
radius in mm	137	147	146	142
<b>Cat. No.</b>	<b>tubes with screw cap</b>			<b>0534<sup>6)</sup></b>



### Adapter

boring Ø x L in mm	26 x 83	29 x 93	38.5 x 92	45.9 x 98
vessels per rotor	4	4	4	4
<b>Cat. No.</b>	<b>4417</b>	<b>4416</b>	<b>1396</b>	<b>0765</b>




2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.  
 3) When using these tubes, bucket 1490 cannot be closed with lid 1492.  
 4) Please remove spacer.  
 5) Tested by TÜV in conformity with DIN EN 61010, section 2-020.  
 6) A rubber stopper for closing the tube for agitating or mixing is available under Cat. No. 0535. The tube may not be centrifuged with the stopper.

## SWING-OUT ROTOR, 4-PLACE | 1324



Rotor	
max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,630
max. capacity	4 x 100 ml
run-up   run-down, braked in sec	27   30
angle   max. noise level	90°   55 dB (A)
<b>Cat. No.</b>	<b>1324</b>

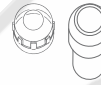

Bucket	
Lid bioseal <sup>5)</sup>	1492
<b>Cat. No.</b>	<b>1490</b>



Vessels															
capacity in ml	0.5	1.5	2.0	1	3	4	5	6	7	9	15	25	50	85	100
Ø x L in mm	10.7x46	11x38	11x38	6x45	10x60	10x88	12x75	12x82	12x100	14x100	17x100	24x100	34x100	38x101	44x100
max. RCF <sup>2)</sup>	2,343	2,415	2,415	2,558	2,594	2,594	2,522	2,522	2,522	2,504	2,504	2,397	2,379	2,576	2,522
radius in mm	131	135	135	143	145	145	141	141	141	140	140	134	133	144	141
<b>Cat. No.</b>	<b>Pediatric</b>	<b>microliter tubes</b>			<b>Rhesus</b>	<b>tubes<sup>2)</sup></b>									

Adapter															
boring Ø x L in mm	11.2x38	11.2x38	11.2x38	6.5x34	10.5x43	10.5x43	13.4x48	13.4x48	13.4x48	17.6x91	17.6x91	25.2x87	35.2x87	38.5x92	45.9x100.5
vessels per rotor	20	20	20	108	36	36	20	20	20	16	16	4	4	4	4
<b>Cat. No.</b>	<b>1351</b>	<b>1351</b>	<b>1351</b>	<b>1339</b>	<b>1343</b>	<b>1343</b>	<b>1383</b>	<b>1383</b>	<b>1383</b>	<b>1329</b>	<b>1329</b>	<b>1330</b>	<b>1331</b>	<b>1396</b>	<b>0761</b>

Vessels														
capacity in ml	1.1-1.4	2.6-3.4	4.9	2.7-3	4.5-5	4-5.5	7.5-8.2	9-10	10	1.6-5	4-7	4-7	8.5-10	15
Ø x L in mm	8 x 66	13x65	13x90	11x66	11x92	15x75	15x92	16x92	15x102	13x75	13x100	16x75	16x100	17x120
max. RCF <sup>2)</sup>	2,540	2,522	2,522	2,522	2,522	2,540	2,540	2,504	2,630	2,522	2,522	2,486	2,486	2,630
radius in mm	142	141	141	141	141	142	142	140	147	141	141	139	139	147
<b>Cat. No.</b>	<b>blood collection / urine vessels</b>													-

Adapter														
boring Ø x L in mm	9x47	13.4x45	13.4x45	13.4x45	13.4x45	15.6x47	15.6x47	17.6x91	17.6x91	13.4x45	13.4x45	16.5x56	16.5x56	17x90
vessels per rotor	28	20	20	20	20	16	16	16	16	20	20	16	16	4
<b>Cat. No.</b>	<b>1457</b>	<b>1383</b>	<b>1383</b>	<b>1383</b>	<b>1383</b>	<b>1459</b>	<b>1459</b>	<b>1329</b>	<b>1329<sup>4)</sup></b>	<b>1383</b>	<b>1383</b>	<b>1348</b>	<b>1348</b>	<b>1347</b>

Vessels												
capacity in ml	15	50	12	25	30	50	10	30	50	85	30	chrome bath tube
Ø x L in mm	17x120	29 x 115	17 x 100	25 x 90	25 x 110	29 x 115	16 x 80	26 x 95	29x107	38x106	44x105	
max. RCF <sup>2)</sup>	2,630	2,630	2,630	2,308	2,630	2,630	2,486	2,415	2,594	2,576	2,504	
radius in mm	147	147	147	129	147	147	139	135	145	144	140	
<b>Cat. No.</b>	<b>tubes with screw cap</b>											<b>0534<sup>6)</sup></b>

Adapter											
boring Ø x L in mm	17x107	30x90	17x80	26x72	26x80	29.5x80	16.5x56	26x83	29x93	38.5x92	45.9x100.5
vessels per rotor	12	4	4	4	4	4	16	4	4	4	4
<b>Cat. No.</b>	<b>1356</b>	<b>1384</b>	<b>6311</b>	<b>1363</b>	<b>1365</b>	<b>6318</b>	<b>1348</b>	<b>4417</b>	<b>4416</b>	<b>1396</b>	<b>0765</b>

## SWING-OUT ROTOR, 4-PLACE | 1324



### Rotor

max. RPM   max. RCF	4,000 min <sup>1</sup>   2,612
max. capacity	4 x 50 ml
run-up   run-down, braked in sec	27   30
angle	90°   56 dB (A)
<b>Cat. No.</b>	<b>1324</b>

### + Bucket

<b>Cat. No.</b>	<b>1398</b>
-----------------	-------------



### Vessels

capacity in ml	5	6	7	9	15	2.6-3.4	2.7-3	4.5-5	4.9	9-10	10	1.6-5	4-7	4-7
Ø x L in mm	12x75	12x82	12x100	14x100	17x100	13x65	11x66	11x92	13x90	16x92	15x102	13x75	13x100	16x75
max. RCF <sup>2)</sup>	2,486	2,486	2,486	2,522	2,522	2,486	2,486	2,486	2,486	2,522	2,522	2,486	2,486	2,397
radius in mm	139	139	139	141	141	139	139	139	139	141	141	139	139	134
<b>Cat. No.</b>	<b>tubes<sup>2)</sup></b>							<b>blood collection / urine vessels</b>						



### Adapter

boring Ø x L in mm	13.4x57.5	13.4x57.5	13.4x57.5	17.5x81	17.5x81	13.4x57.5	13.4x57.5	13.4x57.5	13.4x57.5	17.5x81	17.5x81	13.4x57.5	13.4x57.5	17.5x81
vessels per rotor	20	20	20	16	16	20	20	20	20	16	16	20	20	16
<b>Cat. No.</b>	<b>1486</b>	<b>1486</b>	<b>1486</b>	<b>1482A</b>	<b>1482A</b>	<b>1486</b>	<b>1486</b>	<b>1486</b>	<b>1486</b>	<b>1482A</b>	<b>1482A</b>	<b>1486</b>	<b>1486</b>	<b>1482A</b>

### Vessels

capacity in ml	15	50	12	50
Ø x L in mm	17 x 120	29 x 115	17 x 100	29 x 115
max. RCF <sup>2)</sup>	2,612	2,576	2,522	2,576
radius in mm	146	144	141	144
<b>Cat. No.</b>	<b>tubes with screw cap</b>			



### Adapter

boring Ø x L in mm	17x100	30x98	17.5x81	30x98
vessels per rotor	16	4	16	4
<b>Cat. No.</b>	<b>1483A</b>	<b>1484</b>	<b>1482A</b>	<b>1484<sup>4)</sup></b>

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.  
 4) Please remove spacer.

## SWING-OUT ROTOR, 6-PLACE | 1619




### Rotor

max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,701	
max. capacity	6 x 50 ml	
run-up   run-down, braked in sec	22   25	
angle   max. noise level	90°   54 dB (A)	
<b>Cat. No.</b>	<b>1619</b>	



### Vessels

capacity in ml	15	50
Ø x L in mm	17 x 120	29 x 115
max. RCF <sup>2)</sup>	2,701	2,701
radius in mm	151	151
<b>Cat. No.</b>	<b>tubes with screw cap</b>	

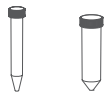

### Adapter

boring Ø x L in mm	17 x 84	30 x 87.5
vessels per rotor	6	6
<b>Cat. No.</b>	<b>1462-A</b>	<b>-</b>



## SWING-OUT ROTOR, 8-PLACE | 1617




### Rotor

max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,469	
max. capacity	8 x 50 ml	
run-up   run-down, braked in sec	22   25	
angle   max. noise level	45°   53 dB (A)	
<b>Cat. No.</b>	<b>1617</b>	



### Vessels

capacity in ml	15	50
Ø x L in mm	17 x 120	29 x 115
max. RCF <sup>2)</sup>	2,469	2,469
radius in mm	138	138
<b>Cat. No.</b>	<b>tubes with screw cap</b>	

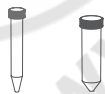

### Adapter

boring Ø x L in mm	17 x 84	30 x 94.5
vessels per rotor	8	8
<b>Cat. No.</b>	<b>1462-A</b>	<b>-</b>



2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.



## SWING-OUT ROTOR, 8-PLACE | 1611



### Rotor

max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,415
max. capacity	8 x 15 ml
run-up   run-down, braked in sec	22   25
angle   max. noise level	90°   53 dB (A)
<b>Cat. No.</b>	<b>1611</b>

### Vessels

capacity in ml	5	6	7	10	15	2.6-3.4	2.7-3	4-5.5	4.5-5	7.5-8.2	1.6-5	4-7	4-7	8.5-10
Ø x L in mm	12/13x75	12x82	12x100	13x100	17x100	13x65	11x66	15x75	11x92	15x92	13x75	13x100	16x75	16x100
max. RCF <sup>2)</sup>	1,914	1,914	2,415	2,415	2,415	1,914	1,914	1,914	2,415	2,415	1,914	2,415	1,914	2,415
radius in mm	107	107	135	135	135	107	107	107	135	135	107	135	107	135
<b>Cat. No.</b>	<b>tubes<sup>2)</sup></b>							<b>blood collection / urine vessels</b>						



### Bucket

boring Ø x L in mm	13x53	13x53	13.2x81	13.2x81	17.5x81	13 x 53	13 x 53	17.5 x 53	13.2 x 81	17.5 x 81	13 x 53	13.2 x 81	17.5 x 53	17.5 x 81
vessels per rotor	8	8	8	8	8	8	8	8	8	8	8	8	8	8
<b>Cat. No.</b>	<b>1131-A</b>	<b>1131-A</b>	<b>1643</b>	<b>1643</b>	<b>1644</b>	<b>1131-A</b>	<b>1131-A</b>	<b>1132-A</b>	<b>1643</b>	<b>1644</b>	<b>1131-A</b>	<b>1643</b>	<b>1132-A</b>	<b>1644</b>

## SWING-OUT ROTOR, 12-PLACE | 1628



### Rotor

max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,683
max. capacity	12 x 15 ml
run-up   run-down, braked in sec	22   25
angle   max. noise level	55° / 60° / 80°   52 dB (A)
<b>Cat. No.</b>	<b>1628</b>

### Vessels

capacity in ml	5	15	2.6-3.4	2.7-3	4-5.5	7.5-8.2	1.6-5	4-7	8.5-10
Ø x L in mm	12/13x75	17x100	13x65	11x66	15x75	15x92	13x75	16x75	16x100
max. RCF <sup>2)</sup>	2,236	2,683	2,236	2,236	2,254	2,683	2,236	2,254	2,683
radius in mm	125	150	125	125	126	150	125	126	150
<b>Cat. No.</b>	<b>tubes<sup>2)</sup></b>		<b>blood collection / urine vessels</b>						



### Bucket

boring Ø x L in mm	13.2x53	17.5x79	13.2x53	13.2x53	17.5x53	17.5x79	13.2x53	17.5x53	17.5x79
vessels per rotor	12	12	12	12	12	12	12	12	12
<b>Cat. No.</b>	<b>1127-A</b>	<b>1621</b>	<b>1127-A</b>	<b>1127-A</b>	<b>1122</b>	<b>1621</b>	<b>1127-A</b>	<b>1122</b>	<b>1621</b>

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.

## ANGLE ROTOR, 8-PLACE | 1418



### Rotor

max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,612
max. capacity	8 x 50 ml
run-up   run-down, braked in sec	36   43
angle   max. noise level	45°   53 dB (A)
<b>Cat. No.</b>	<b>1418</b>



### Vessels

capacity in ml	5	12	1.1-1.4	2.6-3.4	2.7-3	9-10	1.6-5	15	50	15	50	50
Ø x L in mm	12 x 75	17 x 100	8 x 66	13 x 65	11 x 66	16 x 92	13 x 75	17 x 120	29 x 115	17 x 100	29 x 115	29 x 107
max. RCF <sup>2)</sup>	2,182	2,540	2,182	2,182	2,182	2,540	2,182	2,594	2,486	2,540	2,486	2,486
radius in mm	122	142	122	122	122	142	122	145	139	142	139	139
<b>Cat. No.</b>	<b>tubes<sup>2)</sup></b>		<b>blood collection / urine vessels</b>				<b>tubes with screw cap</b>					



### Carrier

boring Ø x L in mm	17.4 x 91	17.4 x 91	17.4 x 91	17.4 x 91	17.4 x 91	17.4 x 91	17.4 x 91	17.4 x 91	30.2 x 92	17.4 x 91	30.2 x 92	30.2 x 92
vessels per rotor	32	32	32	32	32	32	32	32	8	32	8	8
<b>Cat. No.</b>	<b>1467</b>	<b>1467</b>	<b>1467</b>	<b>1467</b>	<b>1467</b>	<b>1467</b>	<b>1467</b>	<b>1467</b>	<b>1468</b>	<b>1467</b>	<b>1468</b>	<b>1468</b>

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.

## ANGLE ROTOR, 6-PLACE | 1620A



### Rotor

max. RPM   max. RCF	6,000 min <sup>-1</sup>   4,226
max. capacity	6 x 85 ml
run-up   run-down, braked in sec	19   22
angle   max. noise level	35°   53 dB (A)
<b>Cat. No.</b>	<b>1620A</b>

### Vessels

capacity in ml	0.5	1.5	2.0	3	15	50	94	7.5 – 8.2	9 – 10	10	8.5 – 10	15	50	50	10
Ø x L in mm	10.7 x 46	11 x 38	11 x 38	10 x 60	17 x 100	34 x 100	38 x 102	15 x 92	16 x 92	15 x 102	16 x 100	17 x 120	29 x 115	29 x 115	16 x 80
max. RCF <sup>2)</sup>	4,105	4,105	4,105	4,105	3,904	4,146	4,226	3,904	3,904	3,904	3,904	3,985	3,985	3,985	3,904
radius in mm	102	102	102	102	97	103	105	97	97	97	97	99	99	99	97
<b>Cat. No.</b>	<b>Pediatric</b>	<b>microliter tubes</b>			<b>tubes<sup>2)</sup></b>			<b>blood collection / urine vessels</b>				<b>tubes with screw cap</b>			



### Adapter

boring Ø x L in mm	11.4 x 39	11.4 x 39	11.4 x 39	11.4 x 39	17.5 x 91.5	35 x 89.3	38.6 x 90.2	17.5 x 91.5	17.5 x 91.5	17.5 x 91.5	17.5 x 91.5	17.5 x 91.5	17 x 106	29.8 x 96.7	-	16.5 x 74
vessels per rotor	24	24	24	24	6	6	6	6	6	6	6	6	3	6	6	12
<b>Cat. No.</b>	<b>1449</b>	<b>1449</b>	<b>1449</b>	<b>1449</b>	<b>1451</b>	<b>1463</b>	<b>-</b>	<b>1451</b>	<b>1451</b>	<b>1451</b>	<b>1451</b>	<b>1451</b>	<b>1466</b>	<b>1454</b>	<b>1646<sup>8)</sup></b>	<b>1448</b>

### Vessels

capacity in ml	30	50	85
Ø x L in mm	26 x 95	29 x 107	38 x 106
max. RCF <sup>2)</sup>	3,824	4,025	4,226
radius in mm	95	100	105
<b>Cat. No.</b>	<b>tubes with screw cap</b>		




### Adapter

boring Ø x L in mm	26 x 85	29 x 92	38.6 x 90.2
vessels per rotor	6	6	6
<b>Cat. No.</b>	<b>1447</b>	<b>1446</b>	<b>-</b>

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.

8) Adapter (Set), 6-place: for conical 50 ml tubes with screw cap.

## — ANGLE ROTOR, 12-PLACE | 1613

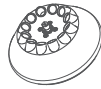

### Rotor

max. RPM   max. RCF	6,000 min <sup>-1</sup>   4,146
max. capacity	12 x 15 ml
run-up   run-down, braked in sec	13   15
angle   max. noise level	35°   55 dB (A)
<b>Cat. No.</b>	<b>1613</b>



### Vessels

capacity in ml	0.5	4	5	6	15	1.1–1.4	2.6–3.4	2.7–3	4.5–5	4.9	7.5–8.2	9–10	10
Ø x L in mm	10.7 x 46	10 x 88	12/13 x 75	12 x 82	17 x 100	8 x 66	13 x 65	11 x 66	11 x 92	13 x 90	15 x 92	16 x 92	15 x 102
max. RCF <sup>2)</sup>	2,777	3,502	3,300	3,300	4,146	3,300	3,300	3,300	4,146	4,146	4,146	4,146	4,146
radius in mm	69	87	82	82	103	82	82	82	103	103	103	103	103
<b>Cat. No.</b>	<b>Pediatric</b>	<b>tubes<sup>2)</sup></b>				<b>blood collection / urine vessels</b>							



### Adapter

boring Ø x L in mm	11 x 35	11.5 x 67.5	13.5 x 60	13.5 x 60	17.7 x 88	13.5 x 60	13.5 x 60	13.5 x 60	17.7 x 88	17.7 x 88	17.7 x 88	17.7 x 88	17.7 x 88
vessels per rotor	12	12	12	12	12	12	12	12	12	12	12	12	12
<b>Cat. No.</b>	<b>2 x 1063-6 (6 pcs.)</b>	<b>6305</b>	<b>1054-A</b>	<b>1054-A</b>	<b>-</b>	<b>1054-A</b>	<b>1054-A</b>	<b>1054-A</b>	<b>-</b>	<b>-</b>	<b>-</b>	<b>-</b>	<b>-</b>

### Vessels

capacity in ml	1.6–5	8	8.5–10	15
Ø x L in mm	13 x 75	16 x 125	16 x 100	17 x 120
max. RCF <sup>2)</sup>	3,300	4,146	4,146	4,146
radius in mm	82	103	103	103
<b>Cat. No.</b>	<b>blood collection / urine vessels</b>			<b>-</b>




### Adapter

boring Ø x L in mm	13.5 x 60	17.7 x 88	17.7 x 88	17.7 x 88
vessels per rotor	12	6	12	6
<b>Cat. No.</b>	<b>1054-A</b>	<b>-</b>	<b>-</b>	<b>-</b>

- 2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.

## — PACKAGES

ROTOFIX 32 A BLOOD TUBE PACKAGE 1*	number of vessels	volume (ml)	size (mm)	RPM	RCF
- 1 x ROTOFIX 32 A centrifuge	1206	20	1.6 - 7	4,000	2,558
- 1 x swing-out rotor, 4-place	1624	16	4 - 10	4,000	2,522
- 4 x bucket	1481				
- 4 x lid (bioseal)	1492				
- 4 x adapter, 5-place	1383				
- 4 x adapter, 4-place	1348				
<b>1206SET1</b>					

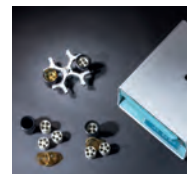

# ROTOFIX 32 A

## Rugged and indispensable

The ruggedly built ROTOFIX 32 A benchtop centrifuge is a medical device used for separating whole blood or blood components of human origin, for example to obtain platelet-rich plasma for autologous therapies. The ROTOFIX 32 A delivers convincing performance thanks to its high capacity and the versatility of its rotor selection.

### — FEATURES

- RPM: 500 - 6,000 min<sup>-1</sup> – adjustable in increments of 100
- Max. RCF: 3,904
- Max. capacity: 8 x 50 ml
- Choice of 9 rotors
- IVDR-conform according to regulation (EU) 2017/746
- Easy operation with keypad
- 2 individual deceleration stages

### — FIELDS OF APPLICATION

- Hospitals
- medical practices
- Dental surgeries
- Orthopedic practices



More information about the control panel can be found on [page 202](#)



\* If you use special PRP kits, please contact us: [info@hettichlab.com](mailto:info@hettichlab.com)



## TECHNICAL DATA

<b>ROTOFIX 32 A</b>	
voltage *)	208 – 240 V 1 ~
frequency	50 – 60 Hz
consumption	300 VA
emission, immunity	EN/IEC 61326-1, class B
max. capacity	8 x 50 ml
max. RPM	6,000 min <sup>-1</sup>
max. RCF	3,904
running time	1–99 min, ∞ continuous run, short cycle mode (impulse button)
dimensions (WxDxH)	366x430x257 mm
weight	approx. 23 kg
noise level	53 dB (A) with rotor 1418
<b>Cat. No.</b>	<b>1207</b>

100 – 127 V 1 ~ / 50 – 60 Hz \*)

1206-01



emission, immunity

FCC class B



\*) Other voltages on request.

## AVAILABLE ROTORS

### SWING-OUT ROTORS

	angle	max. RPM	max. capacity	Cat. No.	page
 swing-out rotor, 4-place	90°	4,000 min <sup>-1</sup>	16x10 ml	<b>1624</b>	68
 swing-out rotor, 4-place	90°	4,000 min <sup>-1</sup>	16x10 ml	<b>1324</b>	68

### ANGLE ROTORS

 angle rotor, 8-place	45°	4,000 min <sup>-1</sup>	8 x 50 ml	<b>1418</b>	69
 angle rotor, 6-place	35°	6,000 min <sup>-1</sup>	6 x 10ml	<b>1620A</b>	69

## SWING-OUT ROTOR, 4-PLACE | 1624



<b>Rotor</b>	
max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,665
max. capacity	16x10 ml
run-up   run-down, braked in sec	22   25
angle   max. noise level	90°   55 dB (A)
<b>Cat. No.</b>	<b>1624</b>



<b>Bucket</b>	
Lid bioseal <sup>5)</sup>	1492
<b>Cat. No.</b>	<b>1481</b>




<b>Vessels</b>	
capacity in ml	4-5.5   4.5-5   4.9   7.5-8.2   7.5-8.2   9-10   10   8.5-10   20
Ø x L in mm	15 x 75   11 x 92   13 x 90   15x92   15x92   16x92   15x102   16x100   31x97
max. RCF <sup>2)</sup>	2,576   2,558   2,558   2,576   2,576   2,540   2,665   2,451   2,308
radius in mm	144   143   143   142   142   142   149   137   137
<b>Cat. No.</b>	<b>tubes / kits</b>

<b>Adapter</b>	
boring Ø x L in mm	15.6x47   13.4x48   13.4x48   15.6x47   17.6x91   17.6x91   17.6x91   17.6x71.5   31x85
vessels per rotor	16   20   20   16   16   16   16   16   4
<b>Cat. No.</b>	<b>1459   1383   1383   1459   1329   1329   1329   1369<sup>4)</sup>   1309</b>

## SWING-OUT ROTOR, 4-PLACE | 1324



<b>Rotor</b>	
max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,630
max. capacity	16x10 ml
run-up   run-down, braked in sec	27   30
angle   max. noise level	90°   55 dB (A)
<b>Cat. No.</b>	<b>1324</b>

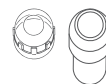

<b>Bucket</b>	
Lid bioseal <sup>5)</sup>	1492
<b>Cat. No.</b>	<b>1490</b>



<b>Vessels</b>	
capacity in ml	1.1-1.4   2.6-3.4   4.9   2.7-3   4.5-5   4-5.5   7.5-8.2   9-10
Ø x L in mm	8 x 66   13x65   13x90   11x66   11x92   15x75   15x92   16x92
max. RCF <sup>2)</sup>	2,540   2,522   2,522   2,522   2,522   2,540   2,540   2,504
radius in mm	142   141   141   141   141   142   142   140
<b>Cat. No.</b>	<b>tubes / kits</b>

<b>Adapter</b>	
boring Ø x L in mm	9x47   13.4x48   13.4x48   13.4x48   13.4x48   15.6x47   15.6x47   17.6x91
vessels per rotor	28   20   20   20   20   16   16   16
<b>Cat. No.</b>	<b>1457   1383   1383   1383   1383   1459   1459   1329</b>

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers.  
 5) Tested by TÜV in conformity with DIN EN 61010, section 2-020.

## ANGLE ROTOR, 8-PLACE | 1418



### Rotor

max. RPM   max. RCF	4,000 min <sup>-1</sup>   2,683
max. capacity	8 x 50 ml
run-up   run-down, braked in sec	36   43
angle   max. noise level	45°   53 dB (A)
<b>Cat. No.</b>	<b>1418</b>



### Vessels

capacity in ml	1.1–1.4	2.6–3.4	2.7–3	9–10	1.6–5	10	50
Ø x L in mm	8 x 66	13 x 65	11 x 66	16 x 92	13 x 75	16 x 100	29 x 107
max. RCF <sup>2)</sup>	2,182	2,182	2,182	2,540	2,182	2,683	2,486
radius in mm	122	122	122	142	122	150	139

Cat. No.	tubes / kits						
	+ 1054-A	+ 1054-A	+ 1054-A	+ 0716	+ 1054-A	+ 0716	+ E2110-A
<b>Carrier</b>							
boring Ø x L in mm	17.4 x 91	17.4 x 91	17.4 x 91	17.4 x 91	17.4 x 91	17.4 x 91	30.2 x 92
vessels per rotor	32	32	32	32	32	32	8
<b>Cat. No.</b>	<b>1467</b>	<b>1467</b>	<b>1467</b>	<b>1467</b>	<b>1467</b>	<b>1467</b>	<b>1468</b>

## ANGLE ROTOR, 6-PLACE | 1620A



### Rotor

max. RPM   max. RCF	6,000 min <sup>-1</sup>   3,904
max. capacity	6 x 10 ml
run-up   run-down, braked in sec	19   22
angle   max. noise level	35°   53 dB (A)
<b>Cat. No.</b>	<b>1620A</b>



### Vessels

capacity in ml	7.5–8.2	9–10	10	8.5–10
Ø x L in mm	15 x 92	16 x 92	15 x 102	16 x 100
max. RCF <sup>2)</sup>	3,904	3,904	3,904	3,904
radius in mm	97	97	97	97

Cat. No.	tubes / kits			
<b>Adapter</b>				
boring Ø x L in mm	17.5 x 92	17.5 x 92	17.5 x 92	17.5 x 92
vessels per rotor	6	6	6	6
<b>Cat. No.</b>	<b>1451</b>	<b>1451</b>	<b>1451</b>	<b>1451</b>

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers.



***LABORATORY EQUIPMENT MAINTENANCE, REPAIR, CALIBRATION AND SALES***

Established in 1987, Henderson Biomedical is the UK's leading laboratory equipment sales and service provider. Our knowledgeable team can provide you with excellent sales advice on a range of different types of laboratory equipment including centrifuges, refrigerators, freezers and heat sealers.

Henderson Biomedical is also able to provide you with first class after-sales service and calibration of your laboratory equipment. We are an **ISO 17025 (UKAS) accredited calibration test laboratory** and our team of Field Service Engineers cover the whole of the United Kingdom.

Please contact us for more information on the types of equipment we supply and the different after-sales services we can offer.

**Henderson Biomedical**

Unit 3, Swan Close  
Croydon CR0 2DZ  
United Kingdom

**Tel: 020 8663 4610**

**For sales enquiries: [sales@henderson-biomedical.co.uk](mailto:sales@henderson-biomedical.co.uk)  
For all other enquiries: [info@henderson-biomedical.co.uk](mailto:info@henderson-biomedical.co.uk)  
[www.henderson-biomedical.co.uk](http://www.henderson-biomedical.co.uk)**

